


Instructions for 'Precision Mount Kit for PD300' (P/N 7Z08220)

Here is a picture of the 'Precision Mount Kit for PD300' (P/N 7Z08220):

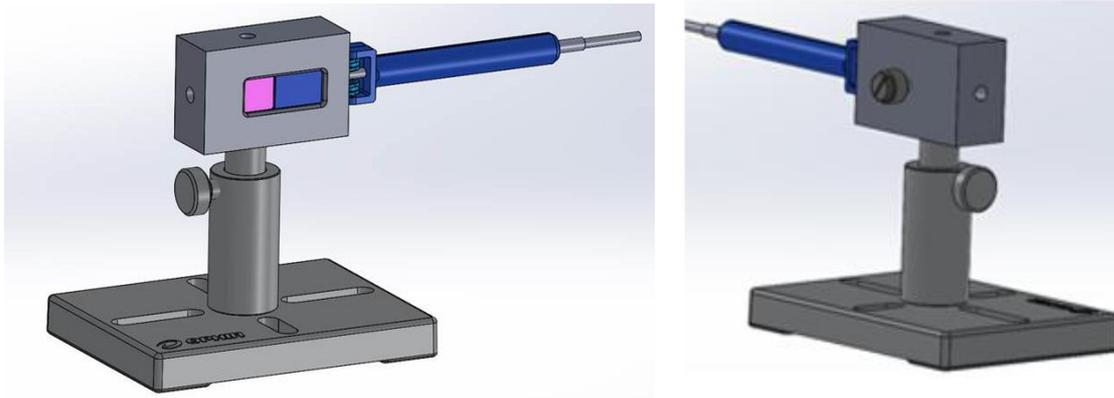

The 'Precision Mount Kit for PD300':


- Assures fixed sensor position in test setup (either vertically or horizontally)
- Allows removal of the sensor (for example re-calibration) and re-insertion to same position

The kit is intended to hold PD300 sensors with an external filter assembled on them. Without the 'external filter' the PD300 sensor will not fit properly into the kit.

Therefore, if the 'external filter' is not needed, the 'filter housing' (see picture below), supplied together with this kit, must be used instead of the 'external filter'. Please make sure to assemble the 'filter housing' on the back of the sensor, so that it will not block the aperture of the PD300 sensor.

Important note: do not assemble the 'external filter' to the back of the sensor, because the screw at the back of the 'Precision Mount Kit for PD300' might damage the 'external filter'.

'Filter housing' (P/N 1G02247) supplied together with this kit:

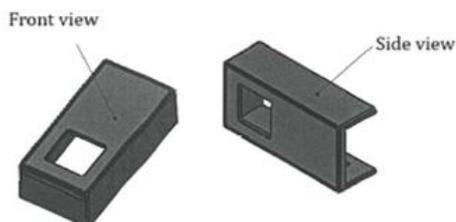

For latest version, please visit our website: www.ophiropt.com